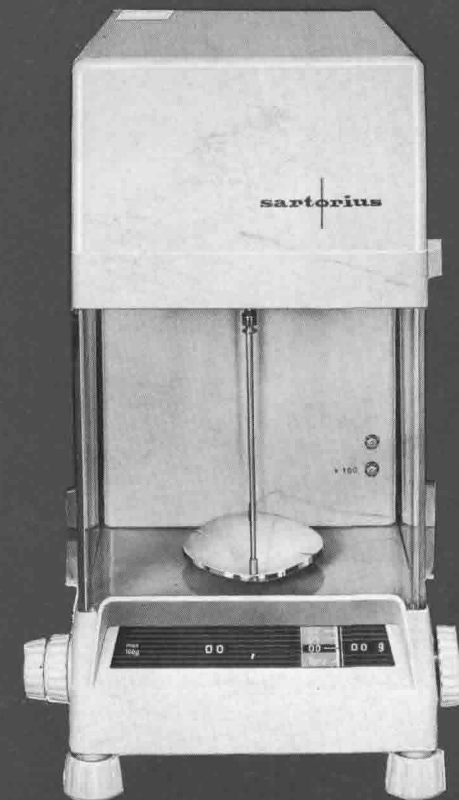
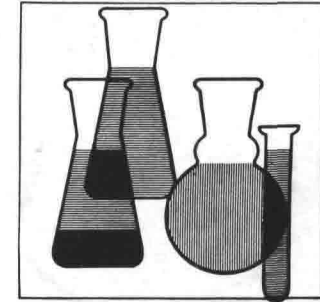
sartorius

Digital Analytical Balances  
Balances Analytiques Digitales  
Balanzas Analiticas Digitales  
Digital Analysenwaagen

# 2400

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After installing the balance we suggest to fix the sticker to the balance or in its neighbourhood.

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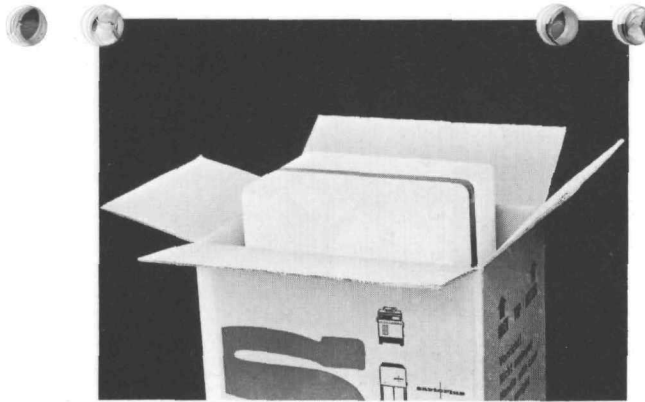
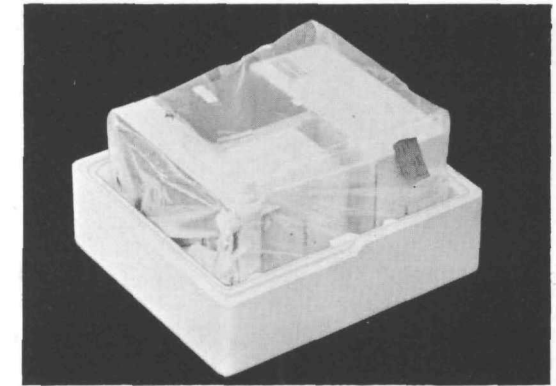
Open Styrofoam insert and take the balance out.

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### Installation

**Note:** Mount balance on a firm floor, sturdy bench or table. Avoid heat radiation and drafts.

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
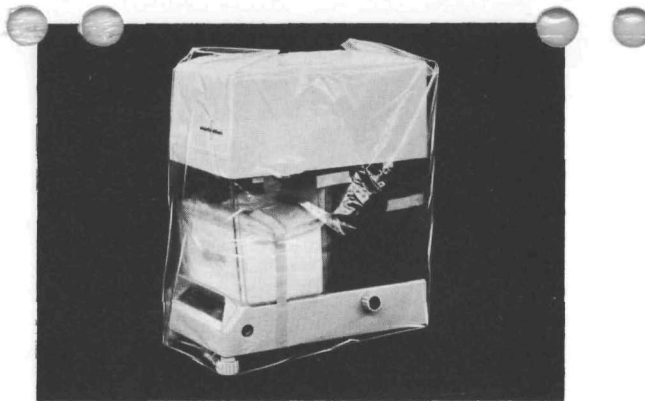
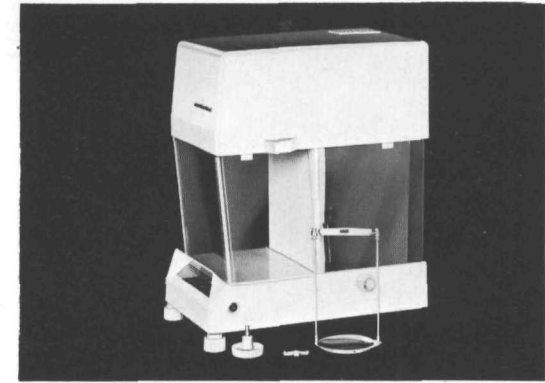
Remove accessories container from the balance. Take out the accessories. Lift hood off balance housing.

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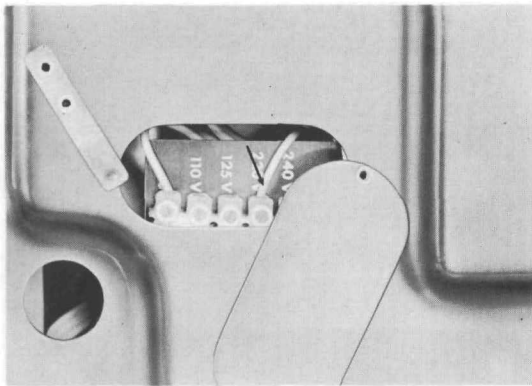
To change the voltage setting:  
 Place the balance with its back on a suitable soft  
 cushion. Open the oblong slider on the bottom  
 cover plate and change connections on the ter-  
 minal strip to the desired voltage.

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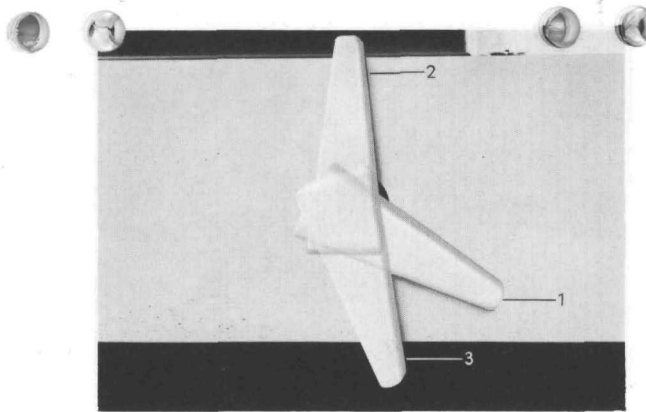Arrestment-Lever on pre-weighing balances

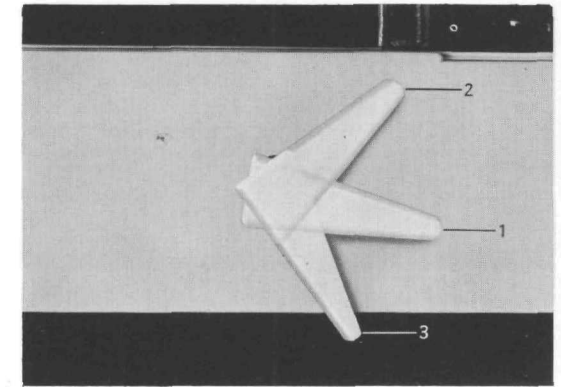
- 1: beam fully arrested
- 2: pre-weighing position
- 3: beam fully released

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Arrestment-Lever on balances without pre-weighing

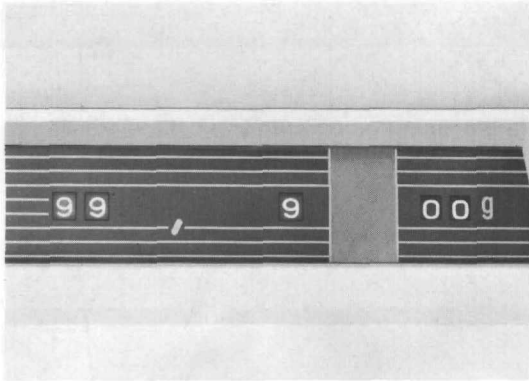
- 1: beam fully arrested
- 2: beam in half release position
- 3: beam fully released

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
With beam-release lever in „arrested“ position.  
Set all weight control knobs to maximum.

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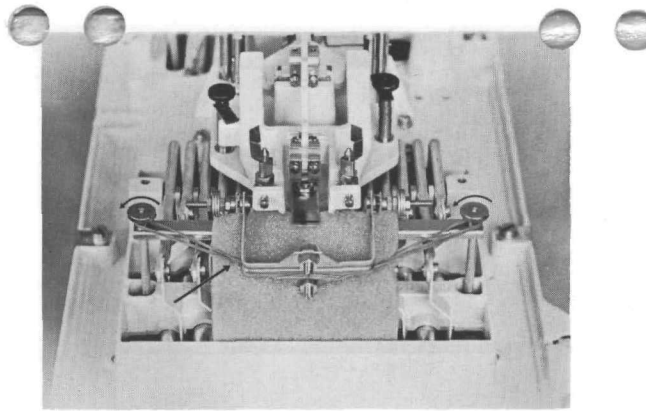
Remove rubber band. Loosen two nuts. Leave  
U-bracket (stirrup) undisturbed.

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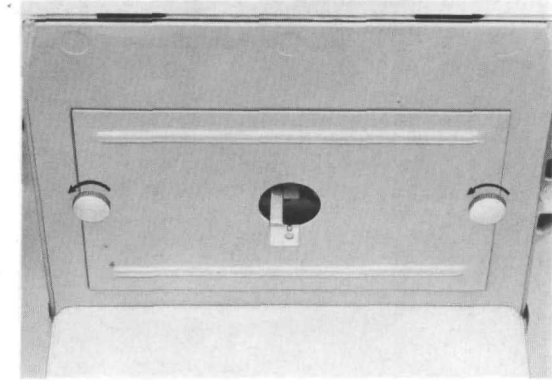
10

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Loosen screws on horizontal dividing panel  
inside the weighing compartment.

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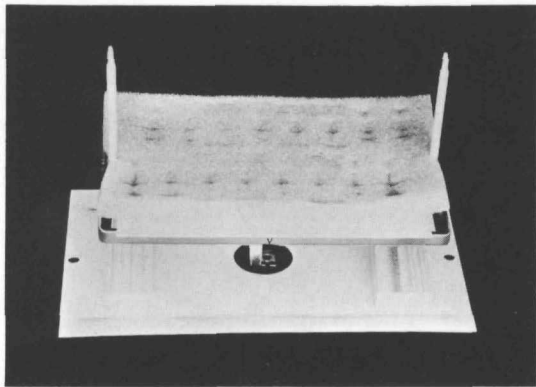
Pull down dividing panel with weight carrier. Remove wooden blocks, cardboard and foam rubber inserts.

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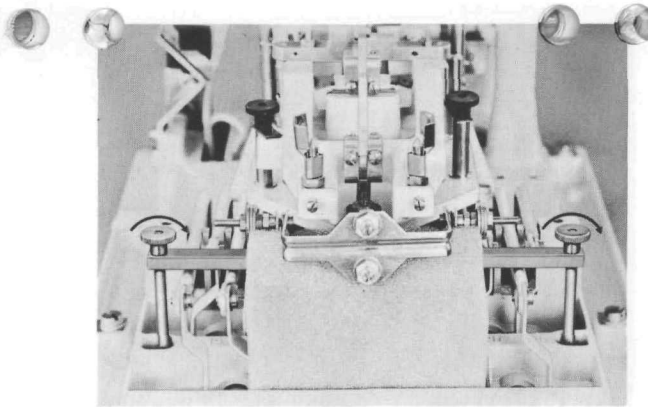
Install weight carrier and attach it to U-bracket. Fasten dividing panel inside the weighing compartment.

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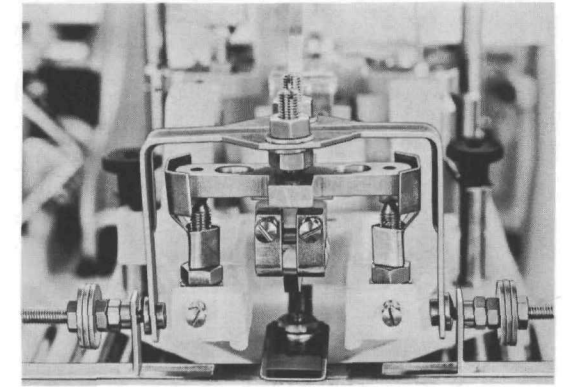
Insert stirrup plate with V-mark facing operator. Remove foam rubber from weight carrier and attach same. Attach pan.

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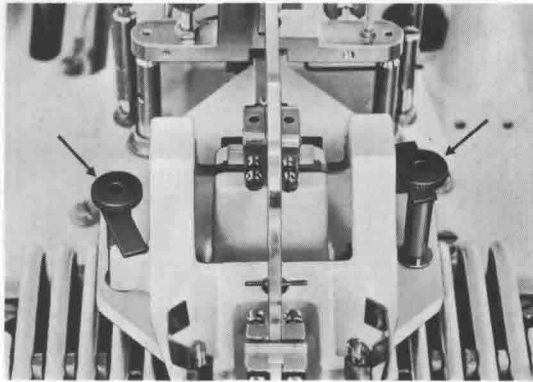
Loosen knurled nuts. Swing leaf springs aside by 90° and re-tighten nuts to prevent accidental displacement of the springs.

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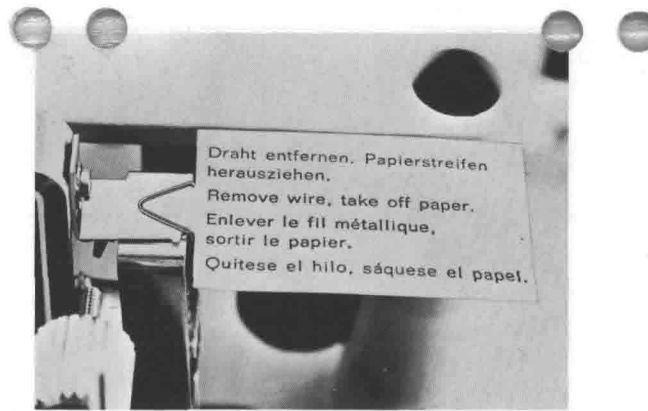
Tare balances:  
Remove clip by pulling out paper strip.

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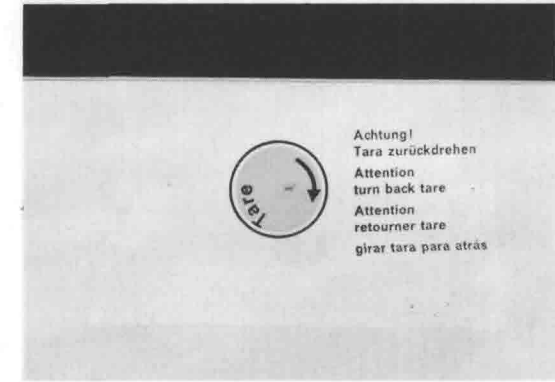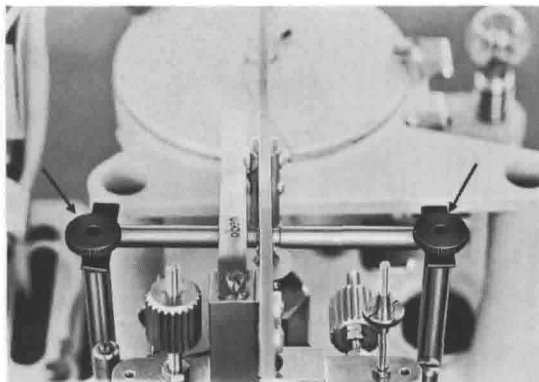
Remove transport arrestment of tare knob. Turn knob counter clockwise to stop.

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
Loosen knurled nuts. Swing leaf spring aside by at least 90°, re-tighten nuts to prevent springs from moving accidentally. Pull out red locking pins and store for possible future use.

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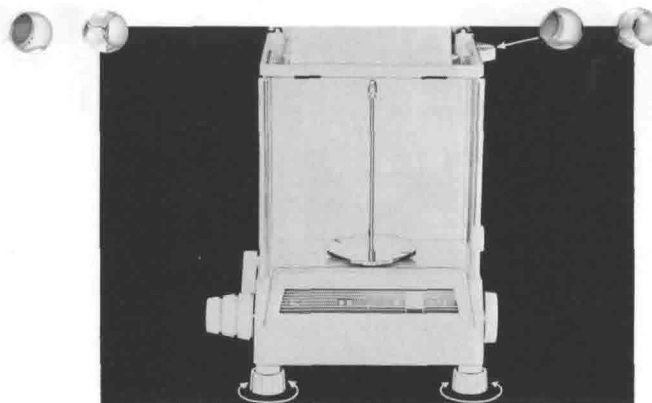
Plug line cord into a.c. outlet. Level balance by adjusting the front feet while observing the spirit level.

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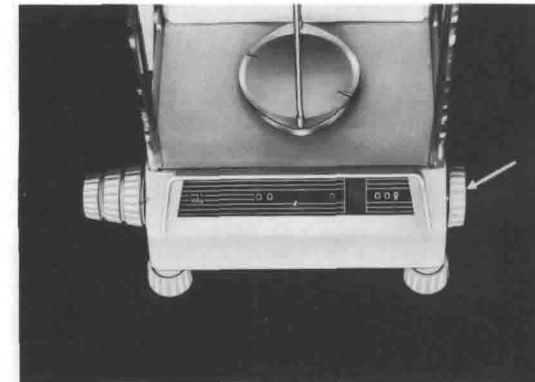
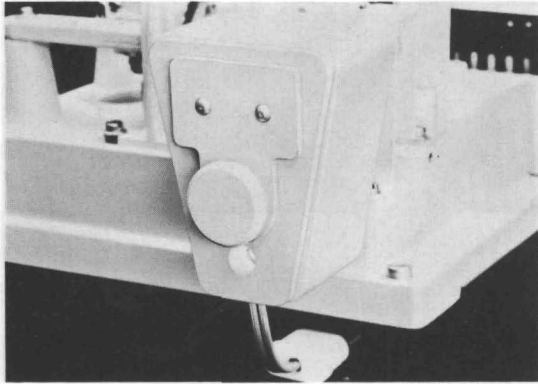
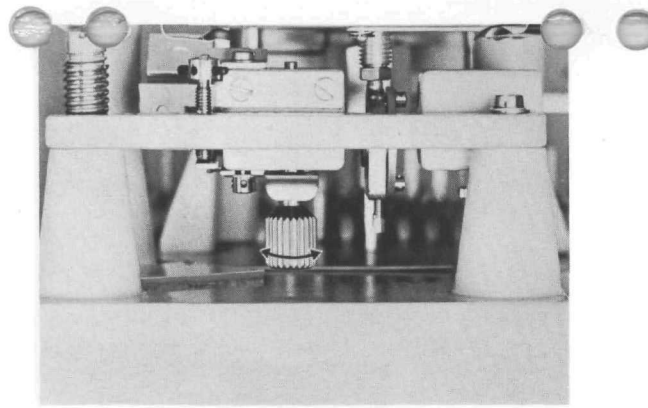
**Checking**  
Release beam arrestment. Focus scale image on projection screen.

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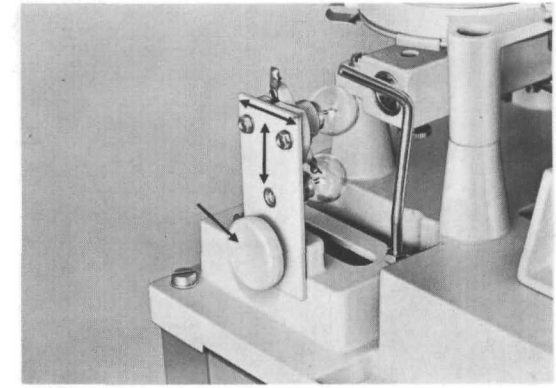
To adjust light intensity. Loosen knurled screw, slide cover plate „sideways“ or „up and down“ until illumination is satisfactory. Tighten screws to fix position of the cover plate; arrest balance.

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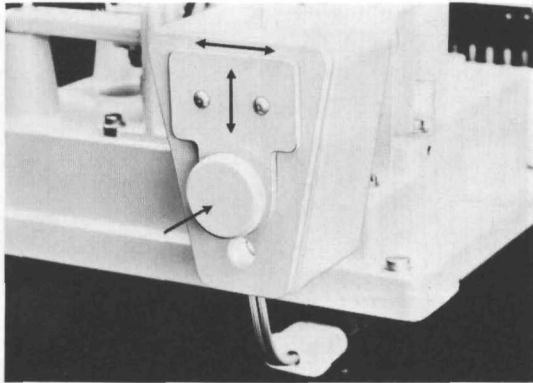
Loosen knurled screw, slide cover plate „sideways“ or „up and down“ until illumination is satisfactory. Re-tighten screws; arrest balance.

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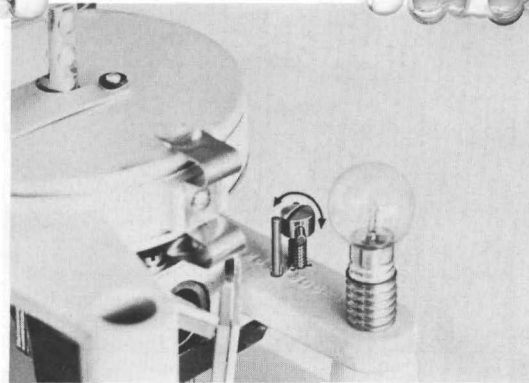
Pre-weighing models: adjust zero point.

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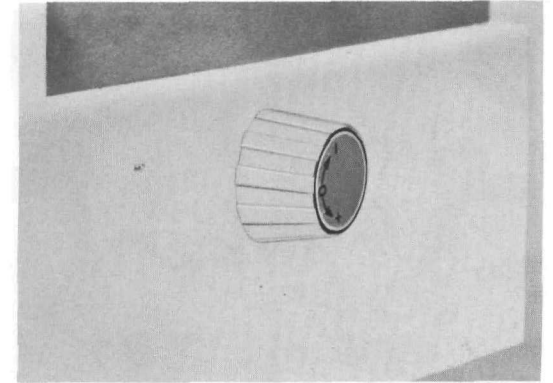
Return hood to housing. Release beam and adjust zero point.

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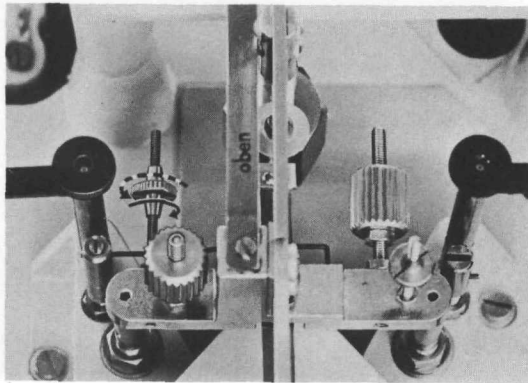1 Umdrehung ca. 10 mg.

If adjusting range is insufficient: Arrest balance. Lift the hood off the balance housing

-  in case of minus deflection
-  in case of plus deflection

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
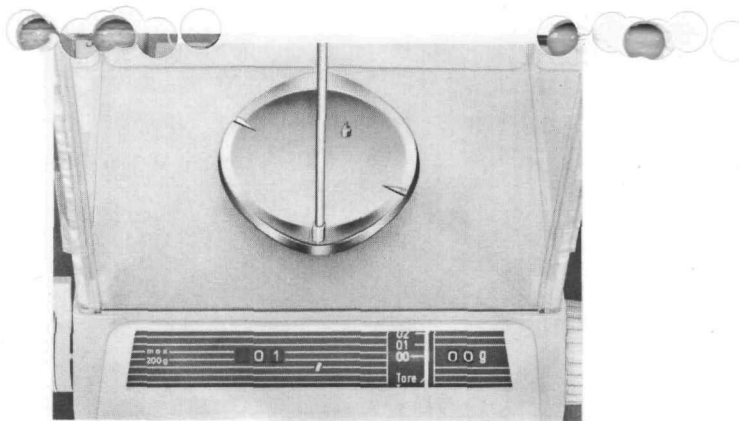
Sensitivity control: Place on the pan a weight which is equivalent to the optical scale range. Switch lowest weight decade to „1“. Release balance.

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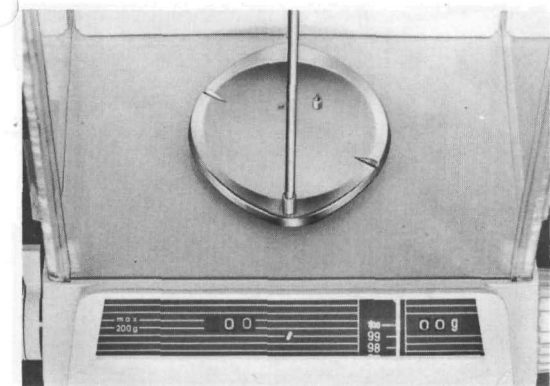
Adjust zero point. Switch weight decades back to „0“. Scale must deflect to „100“.

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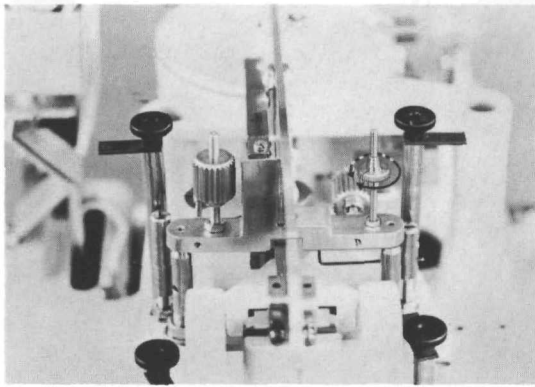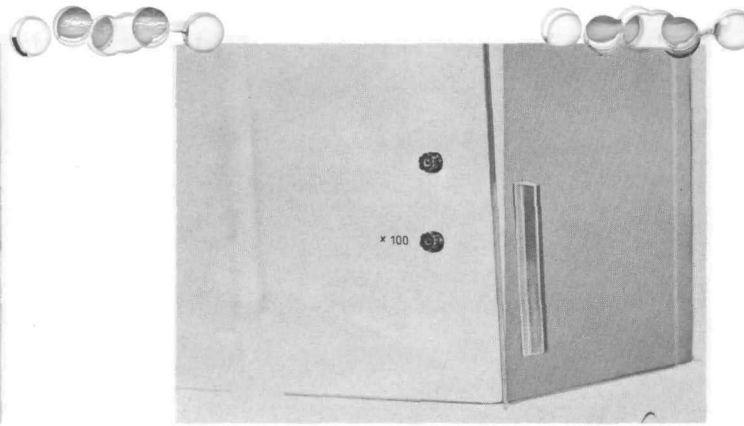
Green Light: Pre-weighing system is operative.

Switch weights to scale value.

Red Light: Beam is in weighing position, do not switch weights now.

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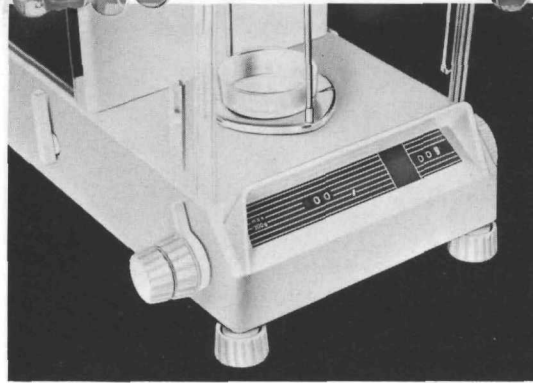
**Weighing, regular**

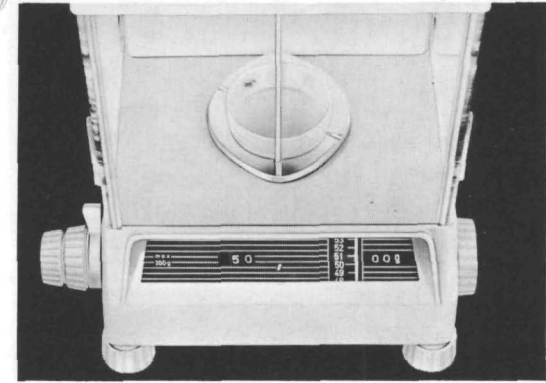
Check zero and adjust setting, if necessary.  
Place load on pan. Move beam release lever UP.

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
Pre-weighing balances:  
Set weight decades in accordance with scale reading.

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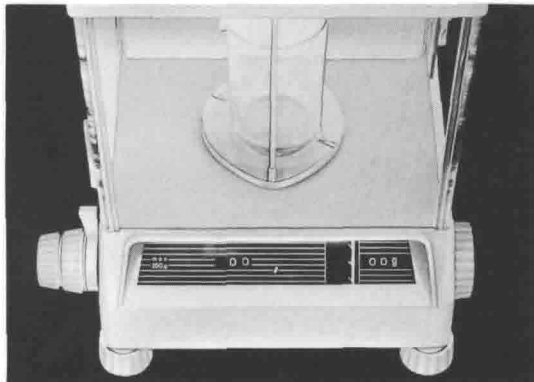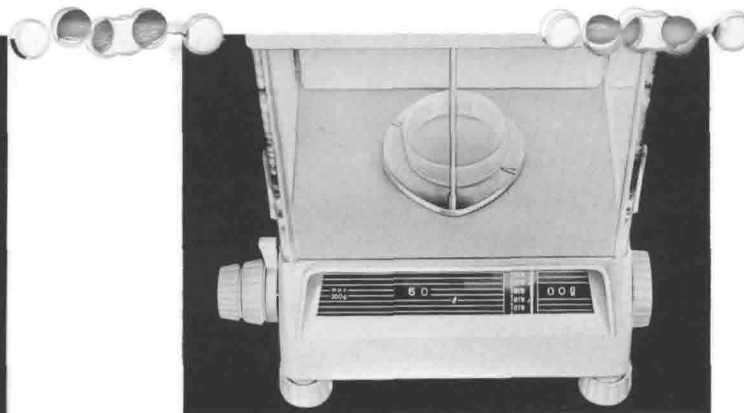
Balances without pre-weighing: Turn knob of the highest decade until the word „Tare“ appears on the screen, then go back one step. Operate the other decades likewise.

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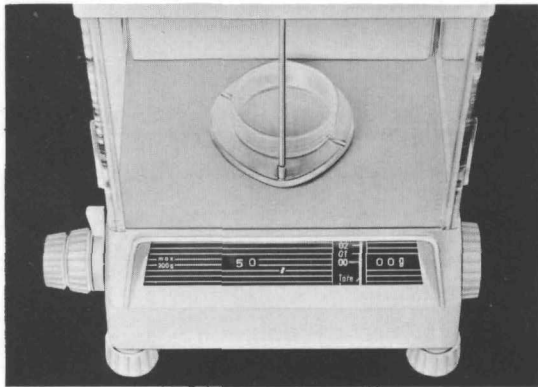
Arrest Beam: After a brief pause, move release lever „down“. When scale comes to a stop, use micrometer next lower division on the scale. Weight indication: 50.3185 g

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
Balances with taring:  
Any reading within the scale range can be tared to  
to „0“.

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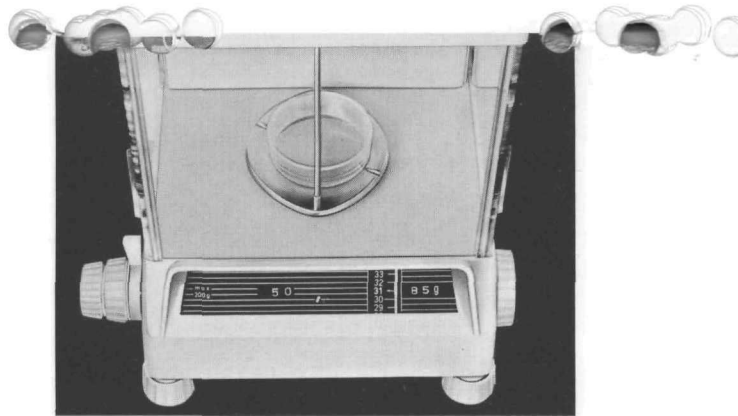**Weighing to pre-set value**

Example:

Tare vessel	50,3185 g
Desired sample weight	<u>2,4850 g</u>
	<u>52,8035 g</u>

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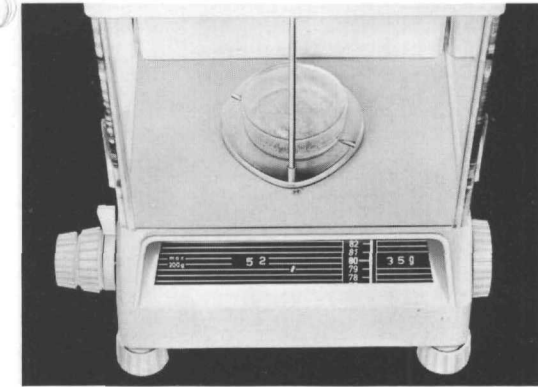
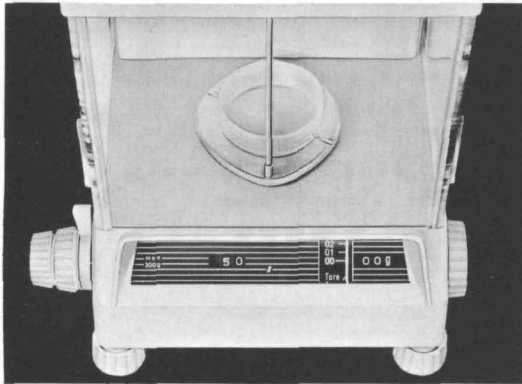
Switch weight decades to „52“.  
Move micrometer to „35“ and add sample until  
scale is at „80“.

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
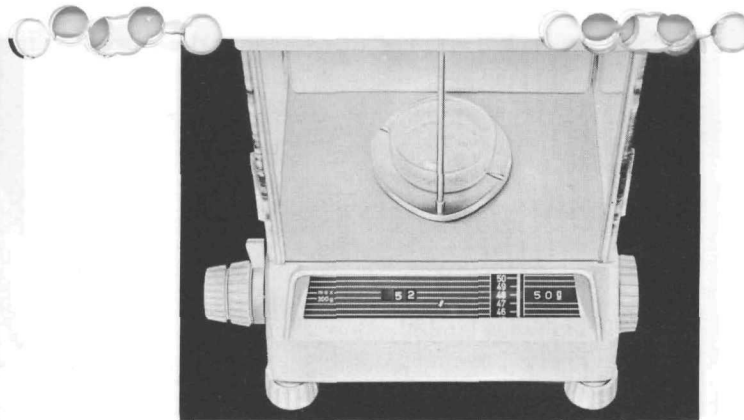
Desired sample weight: 2.485 g  
Switch weight decades to „52“.  
Set micrometer to „50“.  
Add sample material until scale indicates „48“.

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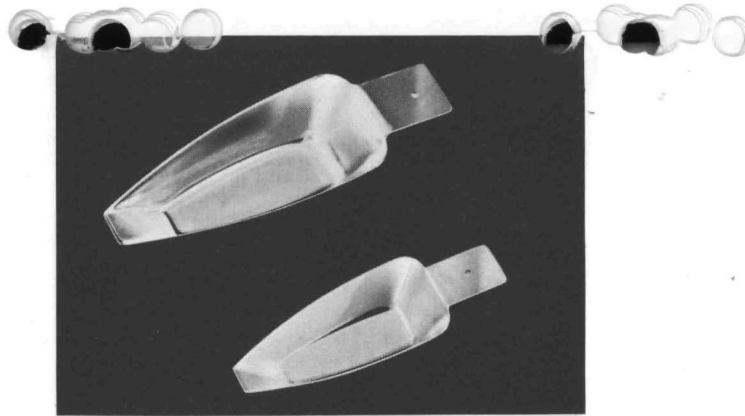Taraausgleichringe 20 g und 50 g.

### Special Accessories

If tared weighing boats are to be used switch weight decades to 20 g or 50 g, respectively.

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
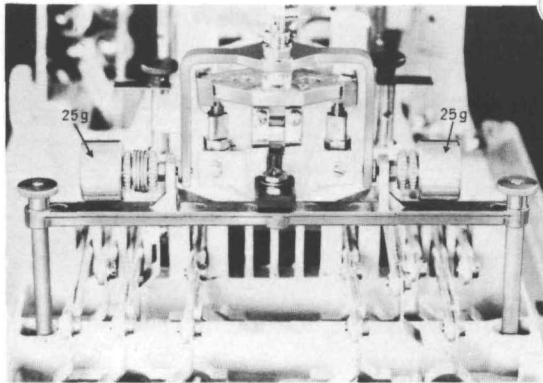
The 50 g tare compensating ring can only be used with balances with two 25 g tare weights.

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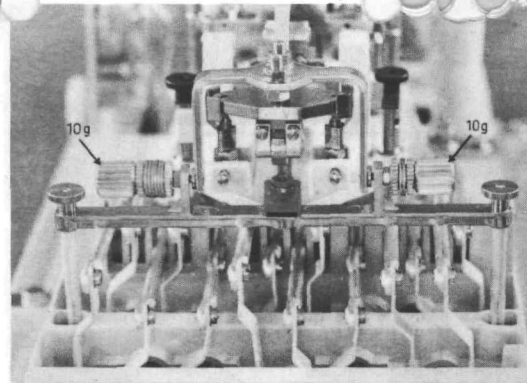
The 20 g compensating ring can be used with all balances equipped with tare weights on the stirrup.

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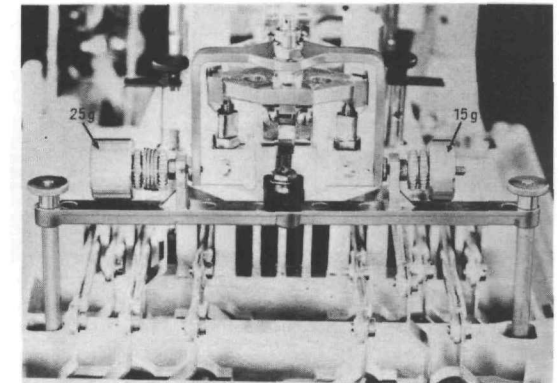When installing the 20 g-tare compensating weights in balances with two 25 g tare weights, exchange these against two 15 g tare weights.

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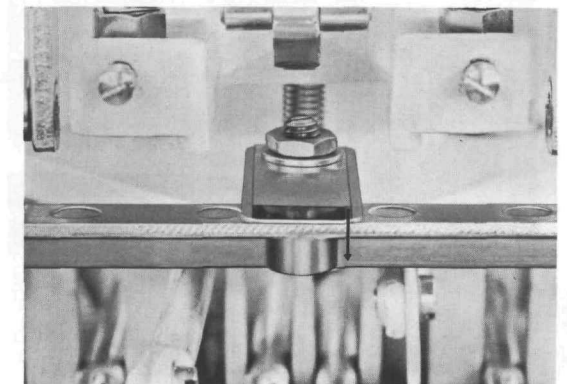
Abb. 53

## Cleaning of the Balance

1. Switch weight decades to maximum settings while balance is in arrested position.

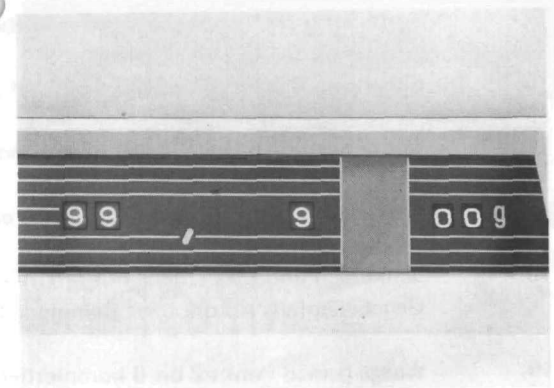
Fig.46

2. Remove pan. Take off stirrup and stirrup plate.
3. Remove the two covers from the damping pot.

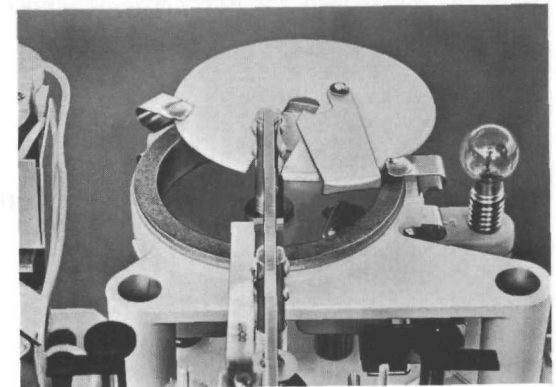
Fig. 47

4. Tare-stirrup: Loosen screw, slightly lift angle bracket.

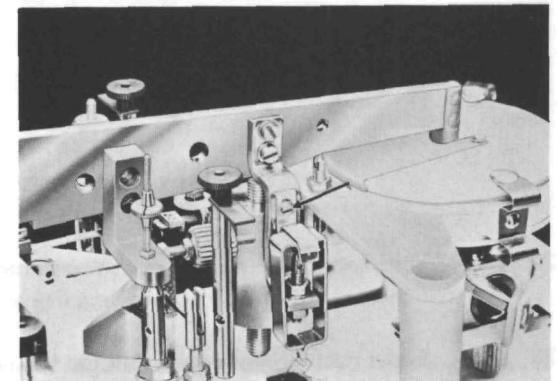
Fig. 48

5. Detach tare hanger and secure properly.

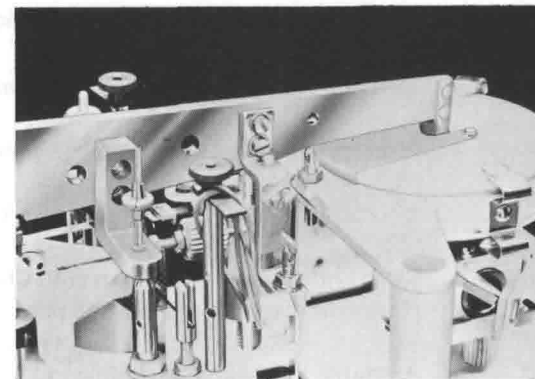
Fig. 49

6. Remove beam.

7. Take weight carrier off as per fig. 10 and 11.

8. Switch one of the decade knobs to „8“. Remove weight from the lowest hook.

9. Switch decade knobs from 7 to 6 to 5 and remove respective weights. Once the weights are removed, return knob to „9“ and proceed accordingly with the remaining decades.

10. Place balance back on a soft cushion. Loosen bottom cover plate. Clean mirror. Reattach bottom cover and set balance on its feet.

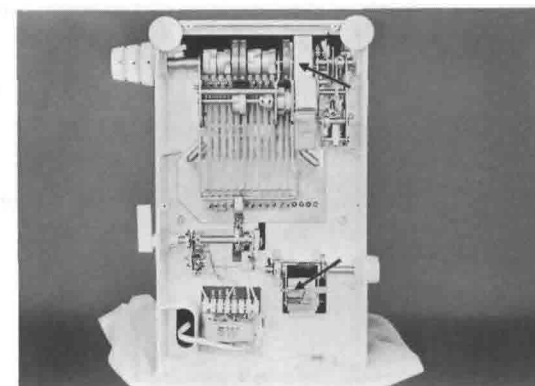
Fig. 50

11. Switch one decade knob to „5“ (others are on „9“).

12. Clean weight with dry Charmois leather and return to proper hook.

13. Weights are marked on one side for identification.

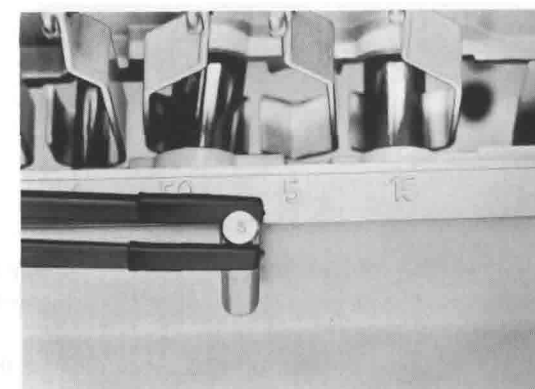
Fig. 51

14. The weight basket is also marked to facilitate proper placing of weights.
15. Switch weight decade from 6 to 7, 8 and 9 and insert the proper weight at each stop.
16. Proceed accordingly with the remaining decades.
17. Insert weight basket according to fig. 12 and 13.
18. Clean knife edges and bearings with Chamois leather and use a magnifier to make sure that all contamination has been removed. If necessary, use alcohol for cleaning.
19. Assemble the balance according to 2 to 6. Adjust zero point according to fig. 25 and 26 check and adjust sensitivity according to fig. 27 - 29.
20. Another check should be performed approximately 2 hours after the various adjustments are completed. A lateral displacement of the scale on the screen not exceeding 2 mm (approximately 1/16") can be corrected by means of this mirror.
21. If the deviation is greater, loosen the optics bracket on one side and move it until the scale is properly positioned on the screen.

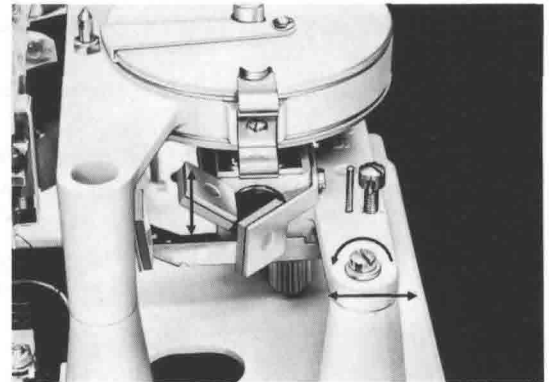
Fig. 52

22. Check pan brake: Push pan all the way back against the panel and release. The pan should stop after 10 to 15 oscillations.
23. Bend the leaf spring down to increase the braking power if oscillation continues longer.

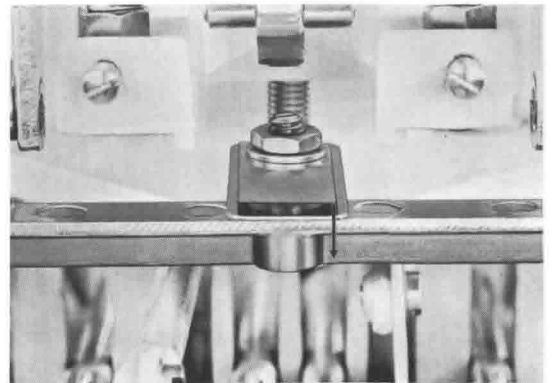
Fig. 53

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AB 2400 · I · 100 · 768 · Mü · Printed in Germany